



## Robust target labeling from small amounts of RNA for Illumina® Expression BeadChip® arrays

The Illumina® Expression BeadChip® arrays are the preferred system for array-based gene expression studies because of their low input requirement, high sensitivity, low cost and high throughput. Two target labeling procedures—the TargetAmp™-Nano and TargetAmp™-Pico Target Labeling Kits for Illumina Expression BeadChip arrays—generate microgram amounts of biotin-antisense RNA (aRNA) from nanogram and picogram amounts of RNA, respectively, and yield high-quality expression analysis results.

Microarray-based transcription profiling studies have provided important advances in our understanding of differential gene expression in heterogeneous cell populations—for example, in cancerous versus normal tissue. The Illumina Expression BeadChips, a popular choice for array-based expression profiling, use biotin-labeled aRNA (sometimes called complementary RNA, or cRNA) as the target. We evaluated BeadChip array results obtained using biotin-aRNA produced by two Epicentre kits: the TargetAmp-Nano and the TargetAmp-Pico Target Labeling Kits for Illumina Expression BeadChip.

### High yields of biotin-aRNA

The TargetAmp-Nano and TargetAmp-Pico kits use an improved Eberwine<sup>1</sup> linear RNA amplification and labeling procedure that generates microgram amounts of biotin-aRNA from nanogram and picogram amounts of total RNA, respectively. The kits have been optimized to maximize the yield of biotin-aRNA and to minimize reaction time and the number of pipetting steps. A TargetAmp-Nano Kit reaction is completed in about 6 h, which enables the user to prepare biotin-aRNA target and begin BeadChip hybridization on the same day.


**Table 1** reports the yield of biotin-aRNA produced from a TargetAmp-Nano kit reaction and from a TargetAmp-Pico kit reaction using various amounts of input total RNA. The actual yield of biotin-aRNA produced from a sample is dependent on the poly(A) RNA content of the total RNA sample, the amount of input RNA and the integrity of the RNA. Both the TargetAmp-Nano and TargetAmp-Pico kits produce sufficient biotin-aRNA for sample replicates or for archiving.

Jim Pease

Jim Pease is at Epicentre (an Illumina company), Madison, Wisconsin, USA. Correspondence should be addressed to J.P. (jim.pease@epicentre.com).

### Genes detected

One of the most important metrics used to evaluate a microarray target labeling method is the number of genes detected by the labeled target. We prepared biotin-aRNA from 100 ng of mouse liver and mouse skeletal muscle RNA with the TargetAmp-Nano kit, and from 50 pg and 100 pg of mouse liver and mouse muscle total RNA with the TargetAmp-Pico kit. The labeled targets were hybridized to the Mouse Ref-8 Expression BeadChip array (Illumina). In a separate experiment, we prepared biotin-aRNA from 500 ng each of RNA extracted from human ovary and testicle using the TargetAmp-Nano kit. The labeled targets were hybridized to the HumanHT-12 v4 Expression BeadChip array (Illumina). The data generated were normalized using BeadStudio (Illumina). The number of genes detected ( $P < 0.05$ ) from each sample is shown in **Figure 1**. The large number of genes detected, even with biotin-aRNA produced from as little as 50 pg of RNA, demonstrates the robust target labeling provided by the TargetAmp kits.

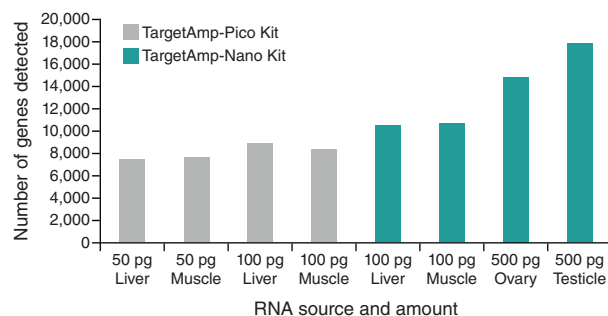

**Figure 1 | Genes detected using labeled target prepared by the TargetAmp-Nano and TargetAmp-Pico kits.** Target RNA from the indicated source was labeled using either the TargetAmp-Nano or TargetAmp Pico kit. The target RNA was hybridized to either the Mouse Ref-8 Expression BeadChip or the HumanHT-12 v4 Expression BeadChip (Illumina).



## APPLICATION NOTES

**Table 1** | Yield of biotin-aRNA from the TargetAmp-Nano and TargetAmp-Pico kits.

Input HeLa cell total RNA	50 µg	100 µg	500 µg	25 ng	100 ng	500 ng
Biotin-aRNA yield (TargetAmp-Nano Kit)				3.1 µg	10.8 µg	75.6 µg
Biotin-aRNA yield (TargetAmp-Pico Kit)	8 µg	19 µg	71 µg			

## Correlation of expression ratios

We also examined the correlation of gene expression ratios using biotin-aRNA produced by the TargetAmp-Nano and TargetAmp-Pico kits. Biotin-aRNA was generated from 100 µg (TargetAmp-Pico kit) and 100 ng (TargetAmp-Nano kit) of RNA extracted from mouse liver and skeletal muscle tissue. There was a strong correlation ( $R = 0.93$ ) between results obtained by the two procedures (Fig. 2), even though there was a 1,000-fold difference in the amount of RNA used for target preparation.


## Conclusions

The TargetAmp-Nano Target Labeling Kit for Illumina Expression BeadChip and the TargetAmp-Pico Target Labeling Kit for Illumina Expression BeadChip kits generate microgram amounts of biotin-aRNA from as little as 50 µg of total cellular RNA. The biotin-aRNA produced detects a large number of genes on the Illumina Expression BeadChip arrays. Expression ratios determined with biotin-aRNA prepared with the TargetAmp-Nano and TargetAmp-Pico kits show high correlation with each other.

## ACKNOWLEDGMENTS

We thank A. Khanna and A. Radek (Epicentre) and I. Lewis and B. Applegate (Illumina, Inc., San Diego, California, USA) for providing data used in this application note.

1. Van Gelder, R.N. *et al.* Amplified RNA synthesis from limited quantities of heterogeneous cDNA. *Proc. Natl. Acad. Sci. USA* **87**, 1663–1667 (1990).

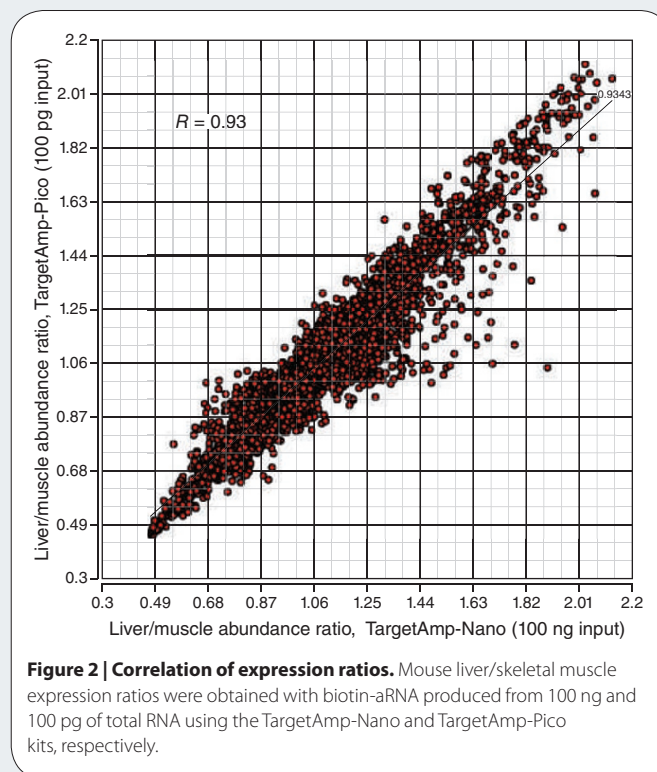

This article was submitted to *Nature Methods* by a commercial organization and has not been peer reviewed. *Nature Methods* takes no responsibility for the accuracy or otherwise of the information provided.